


# Installation of CS2150/CS2350 Bipotentiostat (Ethernet)

## 1. Software installation


- (1) Use the Ethernet cable to connect the computer with the instrument. Power on the instrument.
- (2) Open the “software installation” folder, right click the “Setup.exe”, and run as administrator.

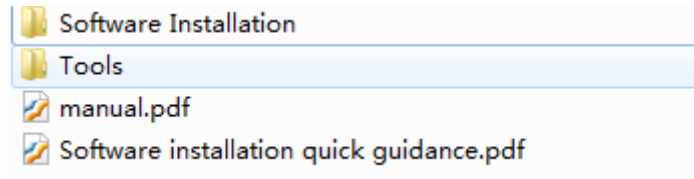

|                 |                 |                |          |
|-----------------|-----------------|----------------|----------|
| CS Studio5      | 2019/8/8 9:41   | 文件夹            |          |
| DotNetFX35      | 2019/8/8 9:41   | 文件夹            |          |
| Drivers         | 2019/8/8 9:41   | 文件夹            |          |
| DriversForXp    | 2019/8/8 9:41   | 文件夹            |          |
| SoConfig        | 2019/8/8 9:41   | 文件夹            |          |
| sxsForWin8      | 2019/8/8 9:41   | 文件夹            |          |
| sxsForWin10     | 2019/8/8 9:41   | 文件夹            |          |
| sxsForWin10-New | 2019/8/8 9:41   | 文件夹            |          |
| clean           | 2017/3/23 9:27  | Windows 批处理... | 1 KB     |
| Setup.exe       | 2019/8/7 15:52  | 应用程序           | 3,769 KB |
| Setup           | 2017/6/29 13:56 | 安装信息           | 1 KB     |
| sxs8            | 2019/5/8 16:57  | Windows 批处理... | 1 KB     |
| sxs10           | 2019/5/8 16:58  | Windows 批处理... | 1 KB     |
| sxs10-New2-x64  | 2019/8/7 13:41  | Windows 批处理... | 1 KB     |
| sxs10-New2-x86  | 2019/8/7 13:42  | Windows 批处理... | 1 KB     |
| sxs10-New-x64   | 2019/5/8 17:58  | Windows 批处理... | 1 KB     |
| sxs10-New-x86   | 2019/5/8 17:58  | Windows 批处理... | 1 KB     |




- (3) As it uses Ethernet communication, no need to install drivers. Firstly, install the .net 3.5. If your computer has already installed it, there will be a prompt as follows.

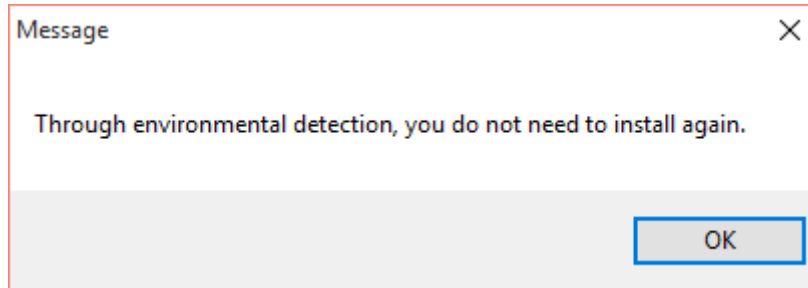


If it has not installed in your computer, just follow the steps to finish installation of .net 3.5.

**Notice:** If your operating system is windows 10 and it appears error prompt when you follow above step, then you have to install .net 3.5 through below tools. Please try one by one till you success. Remember to “run as administrator”.

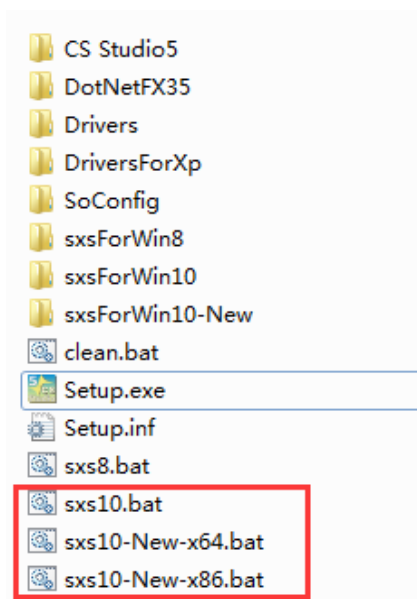


(4) Click install CS Studio5. Created a shortcut is on the desktop.

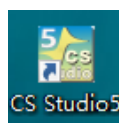


## 2. IP address configuration


Connecting the instrument with computer by Ethernet cable, you should reset the IP address of the computer.

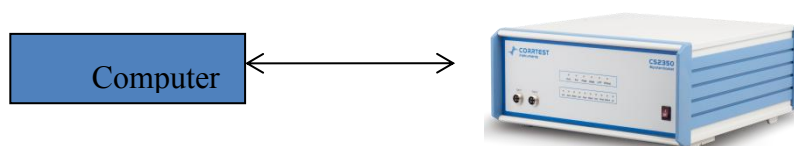

**(1) Find the “Network and sharing center”**

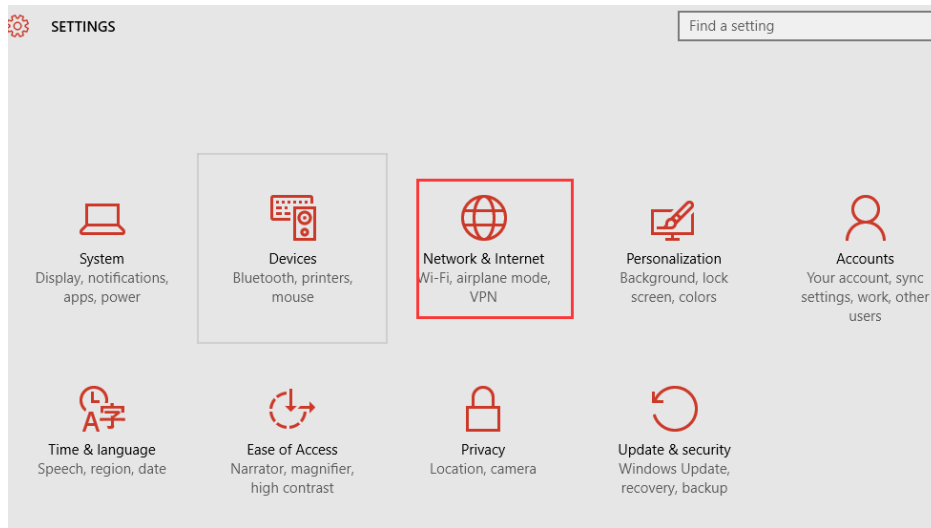


Then, Ethernet → Network and sharing centre

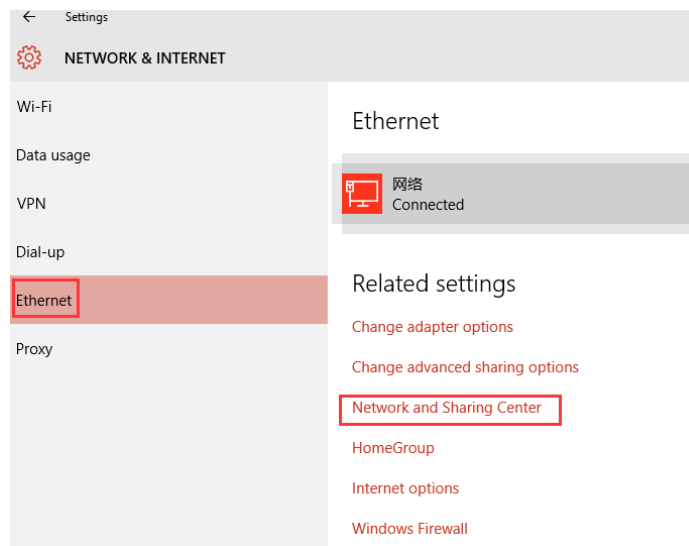



(Or  → Network and sharing center)

**(2) Click “Ethernet” shown as below:**

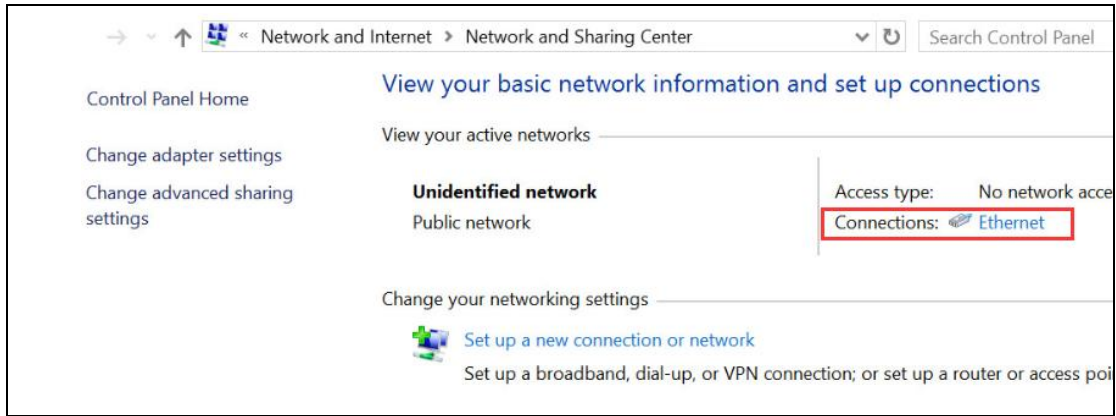


**(3) Click “Properties”, then check “Internet Protocol version 4(TCP/IPv4)”**

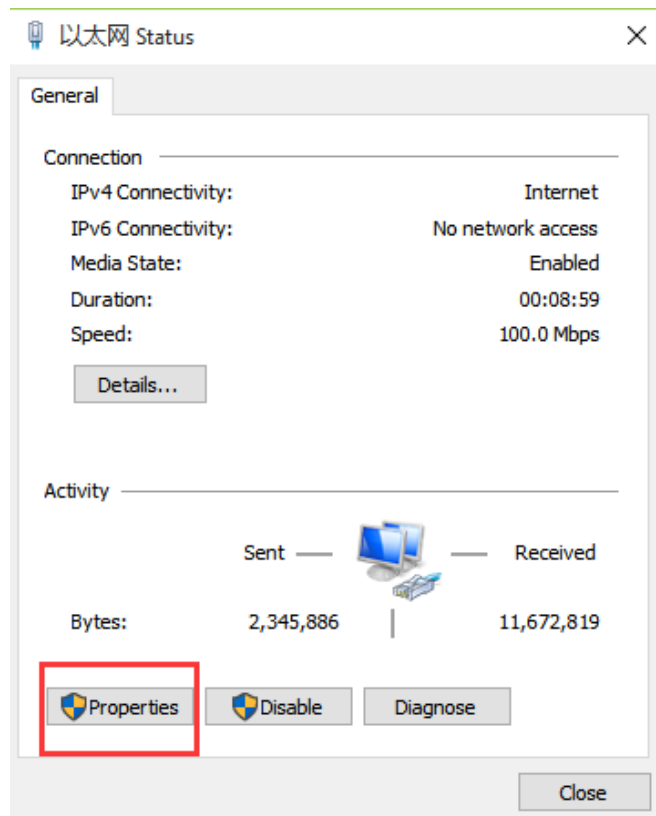


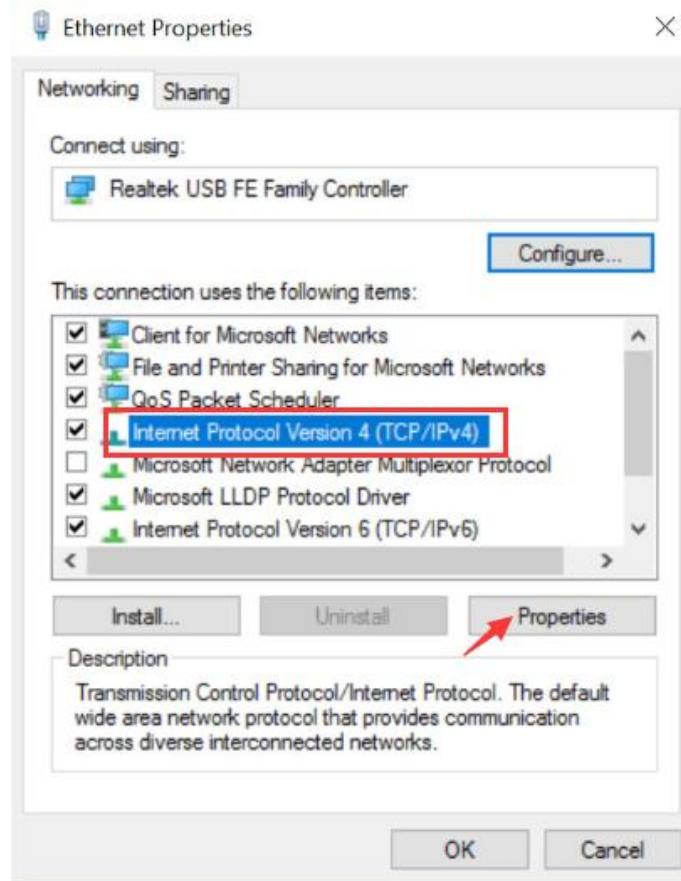

**(4) Fill in the IP address exactly as below:**

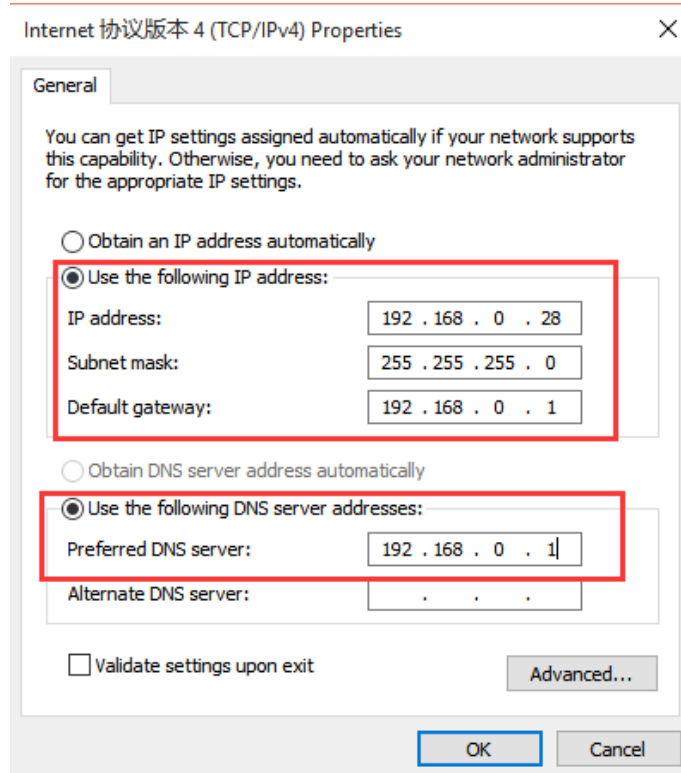

(5) Open CS Studio5 software, “Setup” → Connection → “Ethernet”  
→Manual

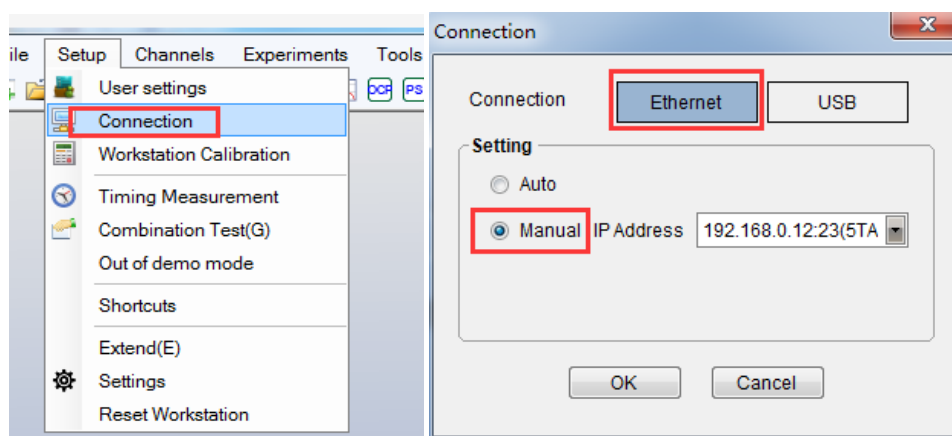

(5) When the bipotentiostat is successfully connected with the computer, there will be model and serial number shown on the software interface, as is shown below (CS2350 for instance).

